

# GamPette™ Precision Powder Pipette

Hand Held Manual Powder Dispensing System

Ideal  
for glove  
boxes



- Repeatedly dispenses powders with a high degree of accuracy.
- Powder is drawn into the tip under vacuum and expelled under light pressure.
- Comprises of a controller unit and handles which are interchangeable and adjustable

Consistent and  
Accurate

GMP Version  
Available

Ergonomic and  
Easy To Use

Typically  
 $\pm 2\%$  RSD

## Precise and Reproducible

Powder dispensing with the **GamPette™** is precise and reproducible. The relative standard deviation achieved on dispensed weights is dependent on the powder properties, however, +/- 2% RSD is typically achieved.

## No Cross Contamination

A rapidly exchangeable, low cost, cellulose filter is incorporated in the delivery line of the pipette to avoid any risk of cross contamination.

## Ergonomic


The lightweight handle allows easy and fatigue free dispensing, ideal for glove box use! The handle is connected to the compact controller unit via a flexible vacuum pipe.

## Versatile

The dispensing volume for each handle is manually adjustable and multiple handle sizes can be used with each **GamPette™** controller unit to cover a wide dispense weight range. The **GamPette™** system is capable of dispensing powder weights from 5 mg to 20 g. Contact us to help you determine the ideal handle size for your target weight range.



## Simple

The pipette is simply operated via a three-position switch (fill - off - dispense) on the hand piece.

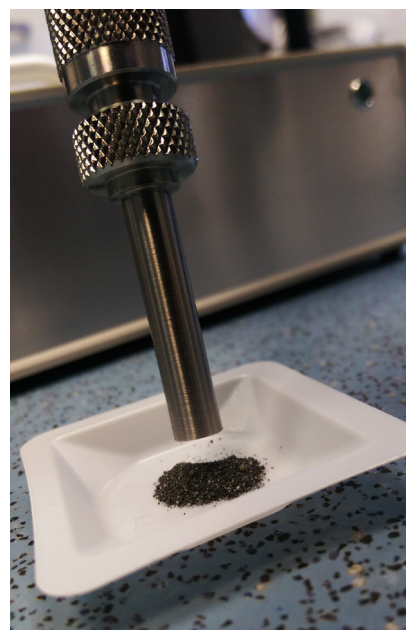

## Typical Gamlen® GamPette™ Dispensing Data

Capable of dispensing powder samples 5 mg to 20 g quickly and accurately.

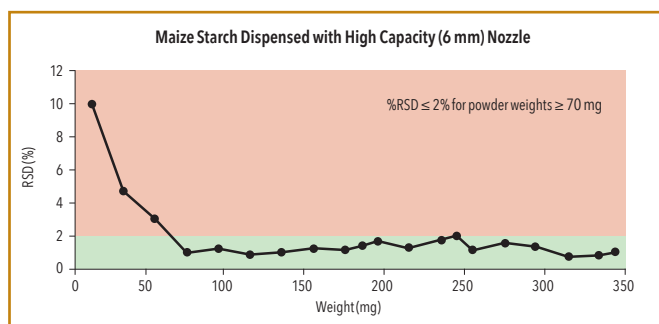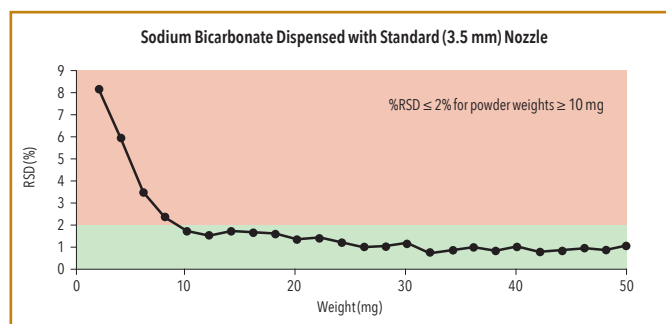

RSD values calculated from 20 measurements

KEY Good Bad

We would be happy to carry out a trial with your chosen material to determine that the expected RSD for your material and target weight range.

## Have a question? Like a quotation? Want to see a demonstration?

Then email [michael.gamlen@gamlentableting.com](mailto:michael.gamlen@gamlentableting.com) or call us now on +44 115 912 4271

Gamlen® Tableting Limited, 38 Croydon Road, Beckenham BR3 4BJ, UK

[www.GamlenTableting.com](http://www.GamlenTableting.com)